
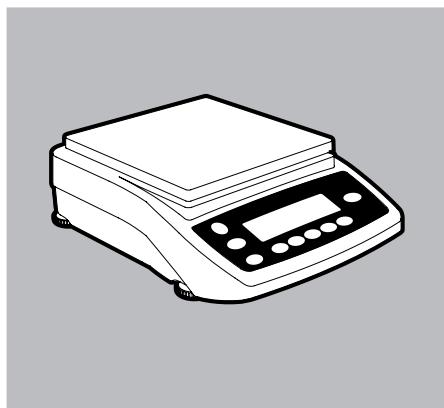
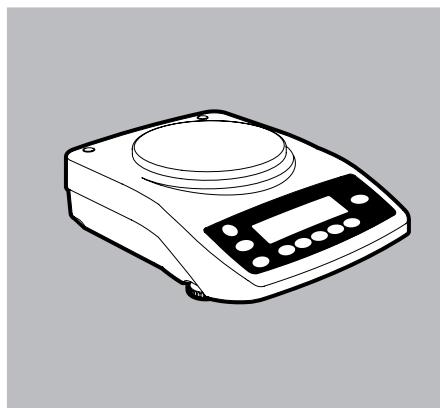
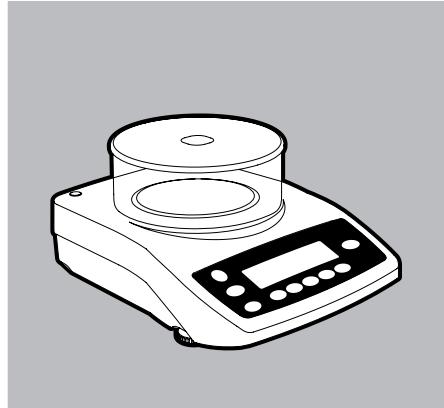
**Operating Instructions | Betriebsanleitung | Mode d'emploi |
Istruzioni per l'uso | Instrucciones de funcionamiento | Handleiding**

Entris

Laboratory Balances | Laborwaagen | Balances de laboratoire
Bilance da laboratorio | Balanzas de laboratorio | Laboratoriumweegschenen

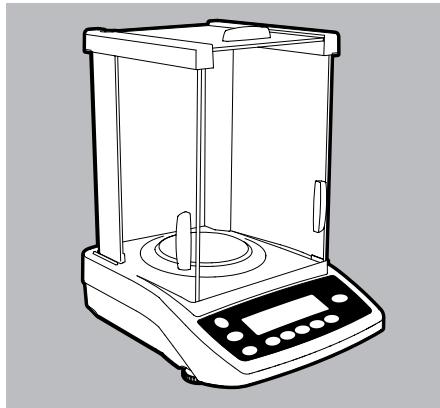98648-019-83

English page 4

In cases involving questions of interpretation, the German-language version shall prevail.

Contents

CD ROM	2
User Information.....	5
Safety Instructions	6
Overview of Display and Operating Elements .	12
Basic Weighing Function.....	13
Calibration and Adjustment.....	14
Configuration (Menu)	17
Declaration of Conformity.....	95

Complete Operating Instructions (Application Programs, etc.):

The enclosed CD ROM contains the instructions as a PDF file for reading and printing.

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Gratis Adobe Reader® van internet te downloaden:
<http://www.adobe.com>

User Information

These notes identify hazards which have a high probability of resulting in death or serious physical injury if not avoided.

These notes identify hazards that can result in moderate or mild injuries if not avoided.

These notes identify hazards associated with the risk of material damage.

Explanation of Symbols

The following symbols are used in these instructions:

- Indicates a required action
- ▷ Describes what happens after you have performed a particular step

Perform steps in the specified order:

1. First action
 2. Second action
 3. ...
- Indicates an item in a list

Conventions for this User Manual:

- The illustrations in these instructions are based on "Standard" balances.

Applications Advice | Technical Support

Addresses for applications advice and technical support can be found on the website at:

<http://www.sartorius.com>

Safety Instructions

Guidelines and General Information

- The balance complies with EU Directives and standards for electrical safety and electromagnetic compatibility*. Improper use or handling can, however, result in damage and/or injury.
Any improper use or operation of the balance, i.e., that is not consistent with the instructions, will result in forfeiture of all claims under the manufacturer's warranty.
- Personnel need to have read and understood these installation instructions, including the safety instructions.
- In the event of use in systems and ambient conditions which have greater safety requirements, you must observe the requirements and provisions applicable in your country.
- Always keep the equipment and balance freely accessible.
Any installation work or balance operation that does not conform to the instructions will result in forfeiture of all claims under the manufacturer's warranty.

* = see "CD-ROM "Specifications" chapter"

Danger of explosion

Do not use this balance in hazardous areas in which explosive materials are present.

Make sure that the voltage rating printed on the AC adaptor is identical to your local mains voltage.

Installation Instructions

Do not operate the balance if the housing or AC adaptor is damaged. Immediately disconnect the damaged device from the power by pulling the plug.

IMPORTANT!

Do not expose the balance, its power supply or accessories supplied by Sartorius to extreme temperatures, aggressive chemical vapors, moisture, shock, vibrations or strong electromagnetic fields.
Observe the conditions of operation described in the "Specifications"!

IMPORTANT!

Note on Installation:
The operator shall be solely responsible for any modifications to the equipment and for connecting any cables or equipment not supplied by Sartorius. Information on operational quality is available upon request from Sartorius. Only use original Sartorius accessories!

Note the IP protection class of the balance and its power supply!
Do not allow liquid penetration.
The protection class specifies the suitability of equipment for various environmental conditions (moisture, foreign bodies).

Before cleaning the AC adapter or the balance: Unplug the power cord.

The balance may only be opened by specialized personnel trained by Sartorius.

Do not open the AC adaptor.

If glass breaks, there is a risk of injury posed by cuts on glass edges.

Lay the cables where they pose no risk of causing someone to trip.

Getting Started

Storage and Shipping Conditions

- Do not expose the balance to extreme temperatures, moisture, shocks, blows, or vibration.

Unpacking the Equipment

- After unpacking the device, check it immediately for any external damage.
- If you detect any damage, proceed as directed on the CD-ROM in the "Care and Maintenance" chapter, "Safety Inspection" section.
- Save all parts of the original packaging for any future transportation. During shipment, please do not leave cables plugged in!

Equipment Supplied

- Balance
- Weighing pan
- Pan support only for models with a round weighing pan
- Mains power supply unit

Additional equipment on the following models:

Entris64-1S, Entris124-1S,
Entris224-1S, Entris64i-1S,
Entris124i-1S, Entris224i-1S

- Sliding panel draft shield
- Shield disk
- Shield plate
- Dust cover

Additional equipment on the following models:

Entris153-1S, Entris323-1S,
Entris423-1S, Entris623-1S,
Entris153i-1S, Entris323i-1S,
Entris423i-1S, Entris623i-1S

- Round glass draft shield
(with shield plate and cover)

Setup


Choose a location that is not subject to the following negative influences:

- Heat (heater or direct sunlight)
- Drafts from open windows and doors
- Extreme vibrations during weighing
- Extreme humidity

Conditioning the Balance

Moisture in the air can condense on cold surfaces whenever the equipment is moved to a substantially warmer place. Allow the device to acclimatize for about 2 hours at room temperature, leaving it unplugged from the power supply.


Installation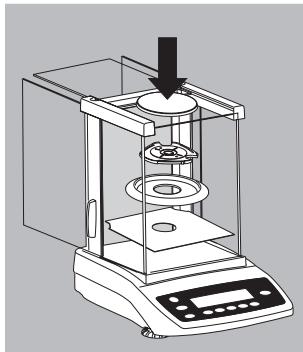

Setting up the Balance


Balances with Sliding Panel Draft Shield

- ▶ Place the components listed below inside the weighing chamber in the order given:
 - Shield plate
 - Shield disk
 - Pan support
 - Weighing pan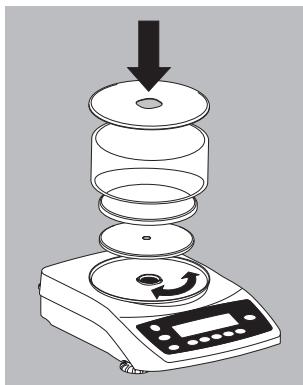

Balances with Round Glass Draft Shield

- ▶ Position the components listed below in the order given:
 - Place lid on top of the balance with the rim facing upwards and rotate until it is securely in place
 - Pan support
 - Weighing pan
 - Glass cover
 - Cover with the rim facing downwards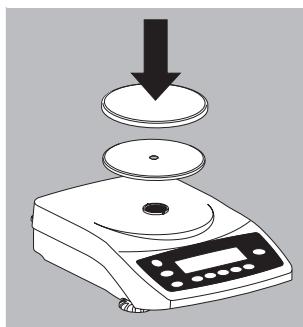


Balances with a Round Weighing Pan

- ▶ Position the components listed below in the order given:
 - Pan support
 - Weighing pan

Balances with a Rectangular Weighing Pan

- ▶ Place the weighing pan on the balance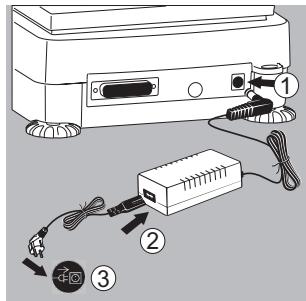

Power Connection | Safety Precautions

- Use only an original AC adapter 6971991
- Insert the right-angle plug into the jack
- Select a country-specific power cable and attach to the AC adapter
- Connect the device to the power

Mains connecting lead:

Item No.	Region Country
6971953	Europe EU (except United Kingdom)
6971954	USA Canada China Japan
6971955	United Kingdom
6971956	Australia, New Zealand
6971957	South Africa
6971964	India
6971998	Brazil
6971999	Argentina
6900931	South Korea


Safety Precautions

Desktop power supply 6971991:

An adaptor rated to Class 2 can be plugged into any wall outlet with no additional safety precautions required.


A ground or earth terminal is connected to the balance housing. The balance housing can be additionally grounded, if required for certain functions.

The data port is also galvanically linked to the balance housing (mass).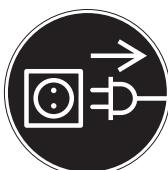


Connecting Electronic Peripheral Devices

- Make sure that the balance is unplugged from the power supply before connecting | disconnecting a peripheral device (printer or PC) to or from the interface port.

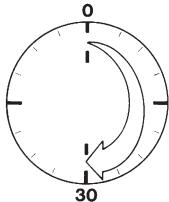Warm-up Time


In order to provide accurate results, the instrument must warm up for 30 minutes. Only after this time will the required operating temperature have been reached.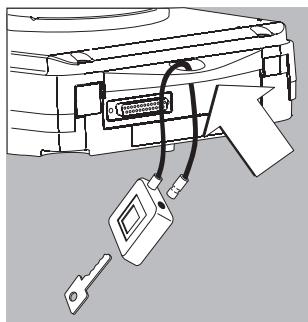

Anti-theft Locking Device

To fasten an anti-theft locking device, use the lug located on the back of the balance.

- Secure the balance at the place of location, e.g., with a chain or a lock.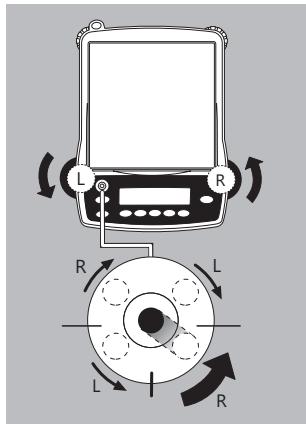

Levelling the Balance

Purpose:

- To compensate for unevenness at the place of installation



Always level the balance again any time after it has been moved to a different location. Only the 2 front feet are used for leveling.

- Screw in both rear support feet (only on models with a rectangular weighing pan).
- Turn the front leveling feet as shown in the illustration until the air bubble is centered within the circle of the level indicator.
- > Normally, several leveling steps are required.
- On models with a rectangular weighing pan:
Screw out both back leveling feet until they touch the setup surface.

Below-Balance Weighing

A port for a below-balance weighing hook is located on the bottom of the balance.

- Not permitted for applications in legal metrology.
- Lift cover plate out of the bottom of the balance.
Attention: Place the balance on its side, do not turn over completely!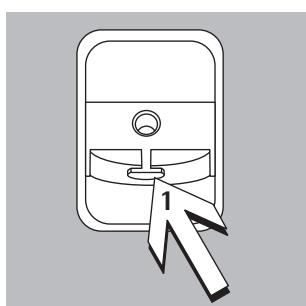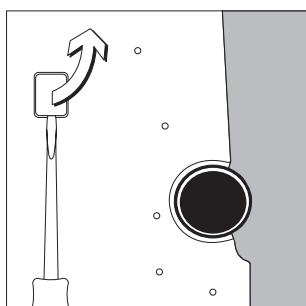


- Secure hook 1: Use a wire, for example, to suspend the sample on the hook.

- Install a draft shield if necessary.

Operation

Overview of Display and Control Panel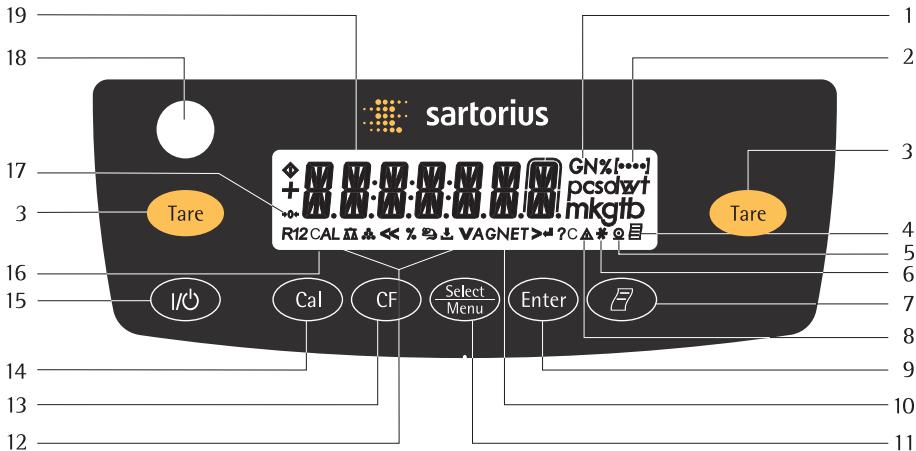

Item	Description
1	Weight units
2	Displays the menu level
3	Taring
4	Symbol for "GLP printing mode active"
5	Symbol for "Printing mode active"
6	Application program active
7	Data output: Press this key to send readout values to the built-in data interface.
8	Calculated-value indicator: not a weight value
9	Start the application program
10	Symbol: Gross or net
11	Select an application program open the operating menu
12	Symbols for active application (ΔΔ, ⚡, %, ⚡, ↓, A, C)

Item	Description
13	Delete (Clear Function) This key is generally used to cancel functions: - Quit application program - Cancel calibration adjustment routine Exit menu
14	Start calibration adjustment routine
15	On Off
16	Symbol: Calibration adjustment function
17	Symbols for "zero range" (verified models only)
18	Level indicator
19	Weight value displayed in selected weight unit Symbol: =<< < < > > <=

Basic Weighing Function

Features

- Tare the balance
- Printing weights

Preparation

- Switch on the balance:
Press the  key
- Tare the balance, if necessary:
Press the  key



► If necessary, change the configuration settings:
see the chapter entitled "Configuration"

► If desired, load the factory settings:
see the chapter entitled "Configuration"

Additional Functions:
► Switching off the balance:
Press 

Example

Simple Weighing

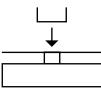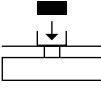Step	Press key	Display Printout
1. Switch on the balance Self-test is performed, followed by automatic initial tare function.		0.0 g
2. Place container on weighing pan (in this example 11.5 g).		+ 11.5 g
3. Tare the balance.		0.0 g
4. Place sample in container (in this example: 132 g).		+ 132.0 g
5. Print weight.		N + 132.0 g

Calibration and Adjustment

Purpose

Calibration is the determination of any difference between the measured value displayed and the true weight (mass) of a sample. Adjustment is the correction of this difference, or its reduction to an allowable level within maximum permissible error limits.

Features

Calibration | adjustment can be performed only when

- there is no load on the balance
- the balance is tared
- the internal signal is stable
- the weight displayed for the sample on the balance must not differ from the nominal weight by more than 2%.

If these conditions are not met, an error message is displayed “ERR 02”.

Following calibration | adjustment, the application program is cleared.

Internal Calibration | Adjustment

Note:

Only for models with the label **Entris...i-1S** !

In the operating menu, select **CAL.JUST. - CAL.INT.** before beginning. The built-in motorized calibration weight located in the housing is applied and removed automatically for internal calibration.

- Select calibration | adjustment:
Press 
- > The internal calibration weight is applied automatically
- > The balance is adjusted | calibrated
- > The built-in calibration weight is removed

Internal Calibration | Adjustment

Note:

Only for models with the label **Entris..i-1S** !

Set the following parameters:

SETUP - BAL.SCAL.- CAL.JUST.- CAL.INT. (Code 1, 1, 9, 4)

The built-in motorized calibration weight located in the balance housing is applied and removed automatically for internal calibration.



Step	Key (or instruction)	Display
1. Tare the balance	(Tare)	0.0 g
2. Start calibration	(Cal)	CAL.INT.
	The internal calibration weight is applied automatically.	CAL.RUN.
3. Adjustment carried out		CAL.END
4. The internal weight is defined automatically		0.0 g

External Calibration

Set the following parameters:

SETUP - BAL.SCAL.- CAL.JUST. - CAL.EXT. (Code 1.1.9. 1)

The required calibration weight is configured at the factory (see CD-ROM "Specifications" chapter).

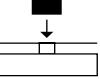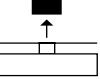Step	Key (or instruction)	Display
1. Tare balance		0.0 g
2. Start adjustment routine		CAL.EXT. 
	Once you store the zero point the display prompts for the required calibration weight (flashing display)	
3. Apply the prompted calibration weight (in this example 5000 g). Weight too low: a minus sign “-” is shown Weight too high: a plus sign “+” is shown		5000.0 g
	The display stops flashing as soon as the weight value is within the defined limit.	
4. Calibration/adjustment executed; then the calibration weight is displayed		CAL.END 
5. Remove the calibration weight		0.0 g

Configuration (Operating Menu)

You can configure the balance; i.e., adapt it to individual requirements.

Functions of the Keys during Configuration:

Symbol	Key	Function
V		Scroll through menu items
>		One menu level lower (use right cursor to scroll through up to 4 menu levels)
↳	  (Press and hold)	Confirm menu item Save settings and exit menu from any position
<<		At the top level: Save settings and exit menu:
<		One menu level higher (left cursor)
[•••]		Indicates menu level

Menu Navigation

Example: Setting the Language

Step	Key (or instruction)	Display
1. Open the menu: Open the menu in weighing mode	 hold	APP LIC.
2. Scroll upward within themenu level; After the last menu code, the first code is displayed again	Repeatedly 	INPUT ... LANGUAG.
3. Select the next menu level (scrolls to the right)	Repeatedly 	ENGLISH °
5. Change setting: Scroll until the desired setting is shown		GERMAN
6. Confirm the menu code; “o” indicates the active setting		GERMAN °
7. Return to the next higher menulevel ▷ Set other menu items as desired	  , 	LANGUAG.
8. Save settings and exit menu	Repeatedly: 	 R2 ALBA 4K 50G VACUUM 1:1 AFB
or		
▷ Exit menu without saving changes		
> Restart your application		0.0 g

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in the form as delivered complies with the essential requirements of the following European Directives and meets the applicable requirements of the harmonized European Standards listed below:

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Dr. Reinhard Baumfalk
Vice President R&D

i.v. Klausre

Dr. Dieter Klausre
Head of International Certification Management

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This declaration certifies conformity with the above mentioned EC and EU Directives, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid. The safety information in the associated product documentation must be observed.

Sartorius Lab Instruments GmbH & Co. KG
Weender Landstrasse 94–108
37075 Goettingen, Germany
Telephone +49.551.308.0
Fax +49.551.308-3289

www.sartorius.com

Copyright by Sartorius, Goettingen,
Germany.

No part of this publication may be reprinted
or translated in any form or by any means
without prior written from Sartorius.

All rights reserved by Sartorius in accordance
with copyright law.

The information and figures contained in
these instructions correspond to the version
date specified below. Sartorius reserves the
right to make changes to the technology,
features, specifications, and design of the
equipment without notice.

Date:
December 2013,
Sartorius Lab Instruments GmbH & Co. KG

Printed in the EU on paper bleached without
chlorine. | W
Specifications subject to change without notice.
Publication No.: WEN6001af131201